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文献资料：Correlations in Infrared Spectra of Nanostructures Based on Mixed Oxides
ISSN 1063-7834, Physics of the Solid State, 2015, Vol. 57, No. 12, pp. 2373–2381. . Pleiades Publishing, Ltd., 2015.

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The IR transmission spectra of the sols containing from 0 to 90 mass % SnO_2 were measured on an FSM 1201 Fourier transform infrared (FTIR) spectrometer (OOO Infracpek, Russia) using an MNPVO36 attachment (with a ZnSe prism)

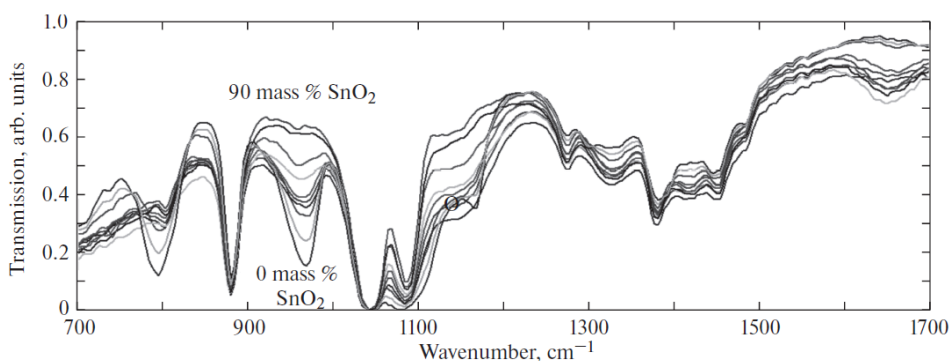
Fig. 1. IR transmission spectra of film-forming sols with different mass fractions of tin dioxide. The SnO_2 content changes from 0 to 90 mass % with a step of 10 mass %.

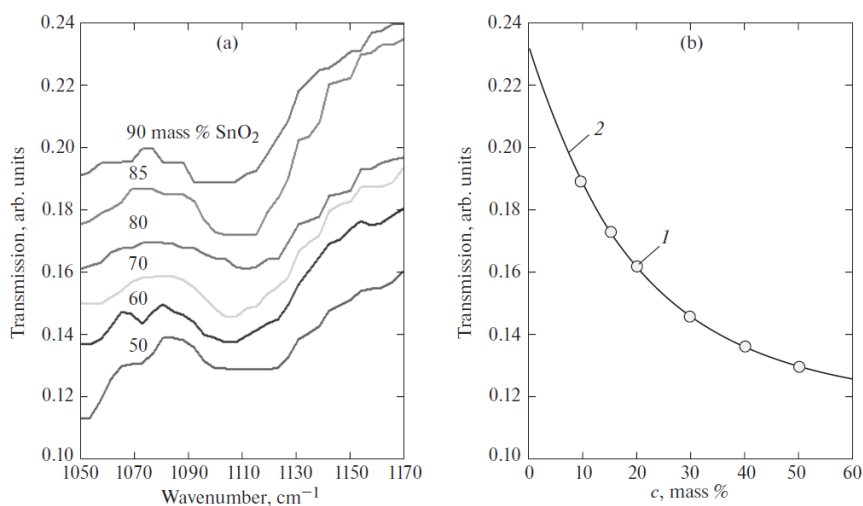
Fig. 6. (a) IR transmission spectra of the nanostructures based on the SiO_2 - SnO_2 oxides. (b) Coefficient of transmission of the infrared light through the studied sols as a function of the SiO_2 mass fraction: (1) experimental data and (2) approximation.

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<p>PZO30 型镜面反射附件</p> <p>镜面反射红外光谱法是一种理想的方法。无需样品预处理就可以进行表面和薄膜的无损检测并测定薄膜厚度。根据镀层厚度选用相应的镜面反射的角度，我供应附件从 10 到 80°。</p>	 A black rectangular device with a silver top panel and a circular opening on the side, used for mirror reflection measurements.
<p>片状糊状样品分析附件</p>	 A collection of laboratory accessories including a mortar and pestle, a small bottle, and several circular sample discs.
<p>液体样品分析附件</p>	 A collection of accessories for liquid samples, including a syringe, a small vial, and several circular sample discs.
<p>气体样品分析附件</p>	 A collection of accessories for gas samples, including a small vial, a syringe, and a circular sample disc.

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